


Maximum Convenience by Advanced Digital Technology

Ultra-Low Temperature Freezer powered by Smart-Lab™ System

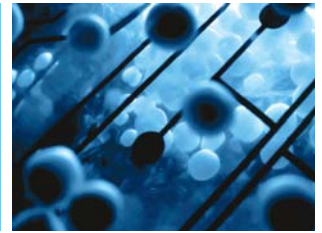

ULT Freezer is evolving like your Smart Phone in hand, and Smart TV at home.

SimpleFreez™ and DuoFreez™ ULT Freezers are the innovation of the freezers. It has been equipped with 7-inch full-touch screen controller and integrated the most extensively applied Android System and Internet Connectivity.

No Button. No Knob. Simply Touch!!

Full Touch Screen Controller and Internet Connectivity!!

The DAIHAN Smart-Lab™ Controller a lot of useful functions like self-diagnostic function, data recording function, and storage management function and so on.

These functions are unique and innovative in the laboratory equipments.

Also this controller is connected with researcher's smart phone, tablet or laptop computer through Internet, by WiFi, so scientists are able to check all the equipments in the laboratory at any time, and at any place.



Smart-Lab™ Controller

- 7 inch full touch screen TFT LCD Controller of "Intuitive Use"
- Data Recording : On-Screen Temperature Chart and USB Storage for Recorded Data
- Internet Connectivity with WiFi
- Remote Monitoring and Control by WiRe™ Service
- Storage Management function : Record the location of samples in the chamber and find easily afterwards, on screen
- Self - Diagnostic Function : Automatic Check & Alarm for the Failures in Key parts

WiRe™ App & Web Service

- Innovative Service by DAIHAN Scientific based on Mobile and Web technology
- Support all different kinds of Web browser by HTML5 Standards
- Available in all the PCs, Laptops and MACs
- Offering Android App and Apple iOS App for Smart phone and Tablet
- Almost same UI as the Touch Screen Controller on the Equipment is available

Smart-Lab™ Controller of SMART Ultra Low Temp. Freezer



- ① USB port : Available Mounted WiFi Module or Data backup and program update of the controller
- ② Display of Present Temperature
- ③ Graph of Recorded Temperature
- ④ Configuration Setting with other functions
- ⑤ Display of Set Temperature
- ⑥ indicating the operation status
- ⑦ Offset Setting
- ⑧ Storage Function to note the sample information in the chamber
- ⑨ Start / Stop
- ⑩ LCO₂ Injection


Features of Smart-Lab™ Controller



Main Display of Controller

» Ergonomic Screen Design

- 7" Full Touch Screen TFT LCD
- Intuitive User Interface
- No-extra Button, Knob or Switch
- All the main Functions and Parameters are displayed on the Main Screen.
- Simple Touch to set the Temperature, Time and other Parameters.

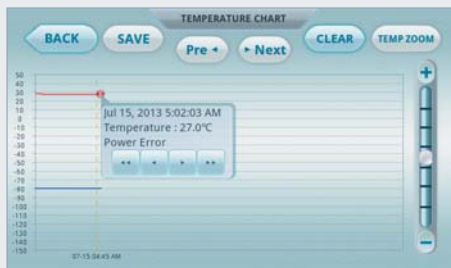

Temperature Chart Display

» Automatic Data Recording

- High Resolution Graphic Display.
- On-Screen Temperature Chart : Check the Temperature Record Graph on the Screen and could see the Details by Zoom-in & Zoom-out enabled at your finger tips.
- The Recorded Data could be transferred to PC by USB Memory.



Error Status Display

» Self Diagnostic Function

- Automatic Check & Alarm for Failures in Key Parts of the equipment.
- Real-time Notification of any Failure in the Equipment to the User(s), to the Service Man and to Supplier, through WiRe™ Service.
- Implemented by LIN and empowered by WiRe™ Service.



Storage Management Function

» Storage Management Function

- Record the location of samples in the chamber and find easily afterwards, on screen.
- Possible to easily and safely manage the location of sample in the chamber.
- Anyone could notice sample's history because it is possible to take memo of sample location.



» Internet Connectivity with WiFi

- Internet Connectivity with WiFi
- WiRe™ App & Web Service : It is possible to monitor and control the freezer, through all smart phone, tablet on Web without restraint in time and place.




» Eco-Friendly Green Freezer

- Dual SHFS™ models offer "Eco-mode", in which the freezer activates the two compressors by turns based on the preset period (8 to 99 days), after reaching down to the set temperature. This specially programmed operation results in dramatically lower energy consumption and much longer lifespan of the compressor.

Maximum Convenience by Advanced Digital Technology

Simple, Efficient and Green - SimpleFreez™ Doubled Security and Power - DuoFreez™

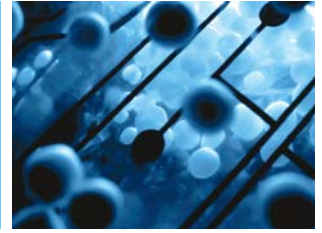

Innovation of Freezing System


DAIHAN SimpleFreez™ System (Simplified Hi-Efficiency Freezing System)

Most of the Ultra-Low Temperature Freezers has two compressors cooperating with each other to drop the internal temperature down to -86°C as shown in the following figure and its overall freezing mechanism is very complicated.


However, our new SimpleFreez™ uses only single compressor. Even though, single set of a compressor, a capillary tube, a condenser and an evaporator is incorporated, it ensures the temperature down to -86°C by an advanced control technology of the compressor and the process of specially mixed refrigerants. SimpleFreez™ overcomes technical complexity of the typical cascaded double compressor system. Simplification of the freezing mechanism and electricity minimizes the noise, power consumption and failure of operation.

High-Efficiency! User Safety! Energy Saving! Long Life time!

Simplified Hi-Efficiency Freezing System Mechanism : Easy Maintenance

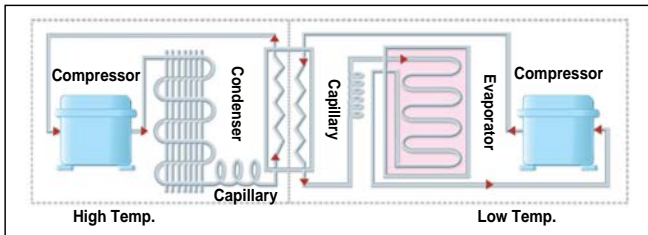


Typical Cascaded Compressor System
Two Freezing cycles contribute Complicated mechanism

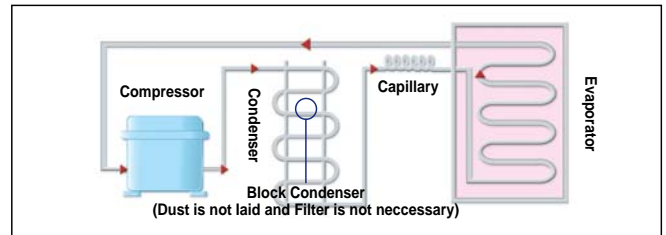

New SimpleFreez™ System
Single freezing cycle works only!!

Doubled Security and Robustness! DuoFreez™

We offer an extraordinary high-stability system powered by two SimpleFreez™ System operating independently. Normally, both of the two compressors are running, and it drops the temperature inside the storage chamber down to -95°C . Even if one SimpleFreez™ System fails in operation by a certain unexpected problem, the other SimpleFreez™ System still works and maintains ultra-low temperature down to -80°C . It means one SimpleFreez™ System works as a 'backup' of the other SimpleFreez™ System. This innovative mechanism realizes doubled stability, robustness and fault tolerance, which are the most important feature of the ULT freezer needed by all the customers.

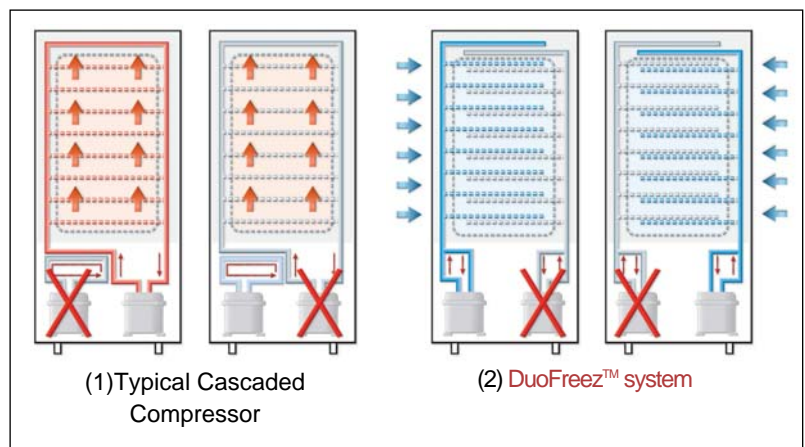

(1) Complicated mechanism

⇒ If one of compressor fails, temperature would go up

(2) Even if one of the two SimpleFreez™ systems fails to run, the other one is still running properly ⇒ Minimum -80°C

▶ The lowest temperature of -95°C is ensured by DuoFreez™ system running independently and simultaneously


▶ Fault Tolerance System

- Realization of doubled stability and robustness

▶ Incomparable Fast and Powerful Freezing

- Fast Recovery of Target Temperature

» Fastest Freezing by DuoFreez™ System

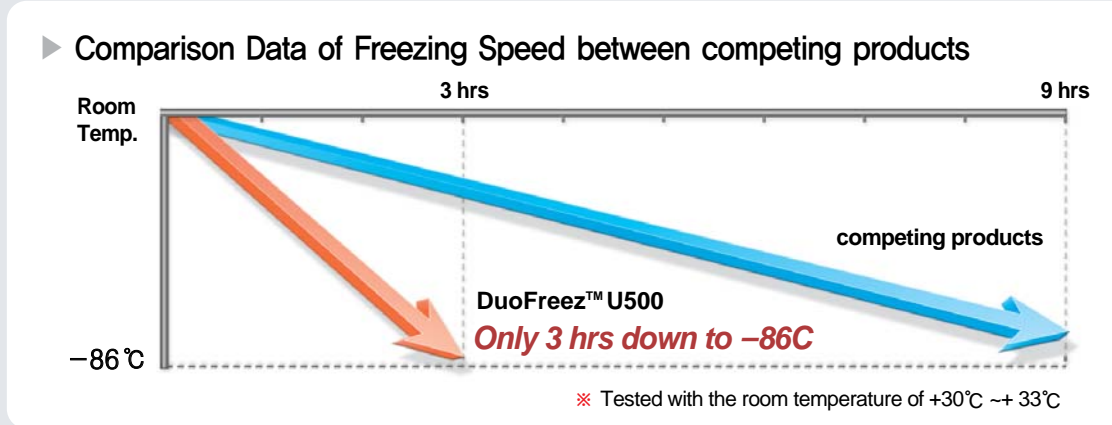

▶ Time Taken to Reach Down to Ultra Low Temperature

Temp.	DuoFreez™			SimpleFreez™				
	U300	U500	U700	U300	U500	U700	WUF-41	WUF-51
	308Lit.	503Lit.	683Lit.	308Lit.	503Lit.	82Lit.	375Lit.	438Lit.
	Upright Type			Chest Type				
- 80 °C				240min.	330min.	150min.	120min.	150min.
- 86 °C	160min.	180min.		270min.	420min.		210min.	270min.
- 90 °C			240min.			270min.		
- 95 °C	270min.	300min.						

※ Tested with the room temperature of 30 ~ 33°C

Real Green Technology for ULT Freezer

» RoHs certified


- The RoHS is the directive on the restriction of the use of certain hazardous substances (Pb, Cd, Hg, Hex-Cr, PBB, PBDE) commonly used in electronic and electronic Equipment (2011/65/EU). All the models of SimpleFreez™ and DuoFreez™ product line are RoHS certified.

» Natural Refrigerants

- The commonly used HFC refrigerants like R-134a, R-404A and R-507, which are potent Greenhouse Gases, will be banned by European law soon. SimpleFreez™ system uses specially mixed refrigerants, which are mostly **Natural Refrigerants**.

» Less Power, Less Money and More Green

- SimpleFreez™ system, extremely simplified freezing mechanism and electricity realize the **lowest power consumption**, leading to **less cost for electricity**. In the end, it also derives **maximum CO2 reduction**, for us to have **more trees** and wider green area.

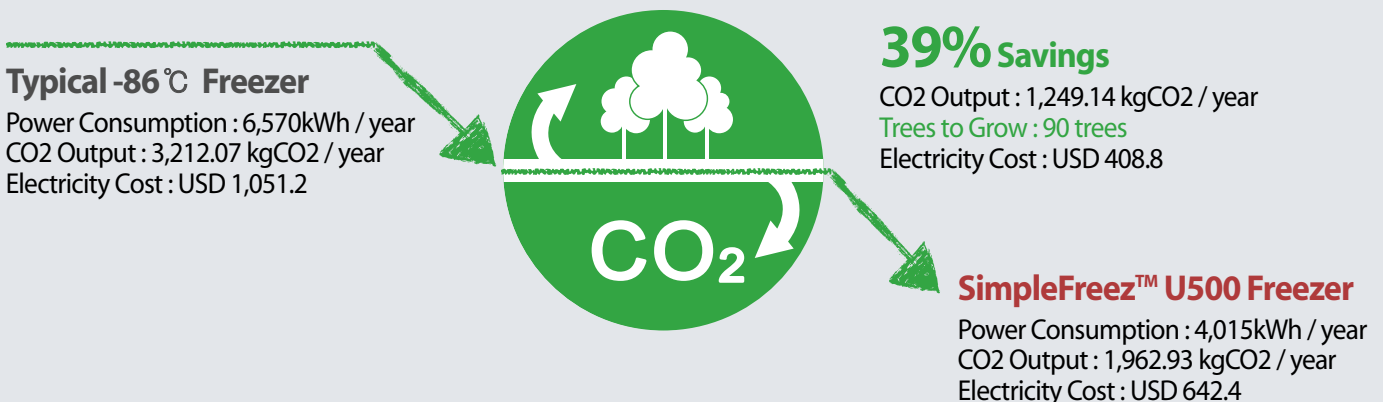

* Calculated by the electricity-specific emission factors sourced from the IEA (2011a).

Ultra-Low Temperature Freezer


Wisdom SMART Ultra



• Robust & Flexible Design of Power Door Handle with Locking function.



• Cable Port for external Temp. sensor

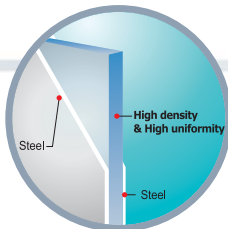

• Insulation Panel packed in housing and door ensures defrosting around outer surface



• Signal Terminal for External Temp. Alarm System

• Filter-Free Mechanism
Significantly Lowers Heat Generated by the Compressor.
Block Condenser (without pins)
Realizes Filter Free Structures.



DuoFreez™ U300



• Automatic Vacuum Dissolver making door opening easy



• Triple for P



Wisdom


DuoFreez™

※ Design of pro




• Temperature Recorder (Optional)
Large and vivid LED

Comparison of Inside formation : SimpleFreez system New technology vs competing products.

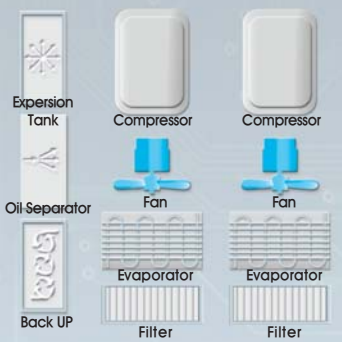

► Complicated Cascaded System

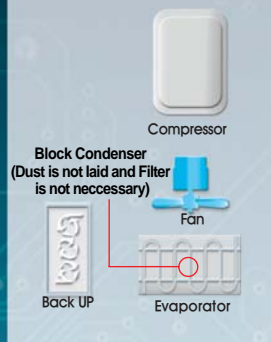

► Innovative SimpleFreez system

- Simplified Hi-Efficiency Freezing System
- Much Simpler
- Oil separator free, Expansion tank free
- Block condenser against typical pin-type condenser - Dust free, Filter free

Low Temp. Freezer



• **Multi-layer Gasket**
- Perfect Insulation.



• **Inner Door Handle**
- easily opened against frosting.



• **Inner door insulation material**




• **Smart-Lab™ Controller**
- Easy setting, Temperature Graph, Storage function (USB interface), etc.

• **CFC-Free**
- HFC refrigerants are used.

• **Maintenance-free cooling device with fully Hermetic High-Capacity compressor.**

• **Alarm**
Built in standard warning and alarm functions.

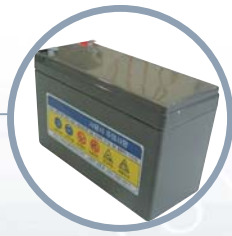


• **Standard Back up Battery System for Data logger and Alarm System.**



C350
duct will be changed

▶ Rack System

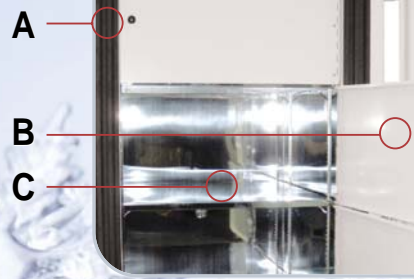

Inside Structure

A High-quality molded thermal insulation material is filled between outer housing and inner chamber, of which structure ensures the stability of the internal temperature and minimizes the frost.

B Four inner doors are equipped for better thermal insulation and efficient use of the storage chamber. They are made of powder coated stainless steel, so it is stronger for anti-frost.

C One piece of internal storage chamber made of high-quality stainless steel is divided into four sub chambers by three shelves, so it's very convenient to efficiently fill the storage boxes inside.



• **Remote Wire/Wireless Alarm System (Optional)**
- Set to be activated when the temperature inside the chamber goes up to -65 °C.
- Main controller signals to RWAS, if connected, at the alarm status, and then RWAS makes phone calls or send messages.
- User could set and store multiple phone numbers and alarm statements.



• **LCO2 back up system (Optional)**
- LCO2 Gas Tank is not included

DAIHAN SimpleFreez™ Personal Digital Ultra-Low Temp. Freezer, Upright Type, -86°C ~ -65°C NEW
 with Smart-Lab™ System (SimpleFreez™ U80 model), Non-Freon Refrigerants, 25-/82-Lit., with Certi. & Traceability



**Personal ULT Freezer "SimpleFreez U80",
with Smart-Lab Controller**

- with Certificate & Traceability : Controlled by Serial Number, Certificate, Delivery-information, and Traceable Data Base System
- Personal type Compact Ultra Low Temperature Freezer
- SimpleFreez™ System can freeze down to -86°C : realize robust and long life time.
- **Smart-Lab™ Controller ("U80" model only)**
- **CE & RoHS Certified**
- **GD(Good Design)-mark ("U80" model only)**
- **PL(Product Liability) Insurance**

- * Personal type Compact Ultra Low Temperature Freezer
- * **SimpleFreez™ System** can freeze down to -86°C : realize robust and long life time.
- * Robust & Flexible design of Door Handle
- * CFC-free Refrigerants are used * External Temp. Sensor Hole
- * Use of Block Condenser removes filter and causes of related failure.
- * Filter-free Mechanism lowers heat generated by the compressor
- * Easy Maintenance and Improved Performance
- * High-quality Insulation Panel and Inner Doors ensure defrosting around outer surface.
- * Built in Standard Warning and Alarm Functions
- * Ergonomic 7" Full Touch Screen TFT LCD (Smart-Lab™ Controller) -only "U80" model
- * WiRe™ App & Web Service - "U80" model only
- * Storage Management Function -only "U80" model
- * Self-Diagnostic Function - "U80" model only
- * Automatic Data Recording - "U80" model only
- * Internet Connectivity with WiFi - "U80" model only
- * RS232 Communication Link and a Remote Control Software are supported - "U80" model only



**Personal ULT Freezer "SimpleFreez U25",
with Digital Feedback Controller**

Specification

Model		SimpleFreez U25	SimpleFreez U80
Capacity		25 Lit.	82 Lit.
Freezing System		SimpleFreez™ System	
Temp.	Range	-86°C ~ -65°C	
	Sensor	at Room Temperature of 30°C and Humidity of 70%	
Compressor		PT100 Sensor	
Condenser Type		Hermetic Type	
Refrigerant		Block Condenser, No Pin, Filter Free Mechanism	
Controller		Natural (CFC Free Refrigerant System)	
Shelves		Digital Feedback Controller	7" Full Touch Screen TFT LCD
Max. Q'ty of Racks		None	Included 1 Stainless steel Shelf
Material	Internal	None, Only 10~15 boxes are applicable	
	External	4pcs of 12(3x4)-Placed Rack (RFU8)	
	Inner Door	Stainless steel (#304)	
Others		Powder Coated steel Plate	
Others		Powder Coated Stainless steel (#304)	
Others		CE & RoHS Certified, GD-mark (SimpleFreez 80) PL(Product Liability) Insurance Touch-Screen TFT LCD Smart-Lab™ Controller (SimpleFreez 80) CFC-Free Refrigerants, RS232 Communication Link Built in Standard Warning and Alarm Functions Remote Control Software (SimpleFreez 80)	
Dimension (w×d×h)	Internal (mm)	200×420×300	340×485×520
	External (mm)	390×627×785	530×687×1,060
Weight		52 kg	60 kg
Packing Size (w×d×h) (mm) & Gross Weight		880×860×1,300	960×760×1,350
Power Consumption		69 kg	102 kg
Power Supply & Cord Plug*		321 Wh	
Power Supply & Cord Plug*		1 Phase, AC 120V, 60Hz or AC 230V, 50 / 60Hz	

* Other Specifications are available upon Customer's Request.

<Continued on next page>

Cat. No	Description	Articles
---------	-------------	----------

<Continued from... **DAIHAN SimpleFreez™** Personal Digital Ultra-Low Temp. Freezer, Upright Type >

Ordering Information


- DH.WUF0032** Upright Personal Type -86°C Freezer, "SimpleFreez U25", -86°C~-65°C, 25 Lit., **230V**
- DH.SWUF0075** Upright Personal Type -86°C Freezer, "SimpleFreez U80", -86°C~-65°C, 82 Lit., **230V**
- DH.WUF0132** Upright Personal Type -86°C Freezer, "SimpleFreez U25", Same as Above, but only **120V**
- DH.SWUF0175** Upright Personal Type -86°C Freezer, "SimpleFreez U80", Same as Above, but only **120V**

Optional Accessories

- DH.RFU8SD** Rack, Sliding Drawer-type, Stainless steel, 12 places, "RFU8SD", w140×d430×h220 mm, for "U80", without boxes
- DH.RFU8SA** (1) Rack, Side Access-type, Stainless steel, 12 places, "RFU8SA", w140×d430×h220 mm, for "U80"
- C11.001** (2) Hard Board Box, with Hard Board Divider, 81 Hole, 127×127×h52 mm
- C11.002** , with PolyPropylene Divider, 81 Hole, 127×127×h52 mm
- DH.KT622** (3) Temp. Chart Recorder, "KT622", with "C657" Chart Paper
- DH.C657** (4) Chart Paper, "C657", 60 Charts, for 420 days recording
- DH.P222** (5) Replacement Pens, "P222", 6ea Red color, for "KT622" Chart Recorder
- DH.LCO2B** (6) LCO2 Back up System, "LCO2B"
- DH.RWAS** (7) Remote Wire/Wireless Alarm System, "RWAS"

Validation (IQ, OQ) Service

- DH.WUF30032** Validation Service(IQ, OQ), "UFV32", for "SimpleFreez U25"
- DH.SWUF30075** , "SUFV75", for "SimpleFreez U80"

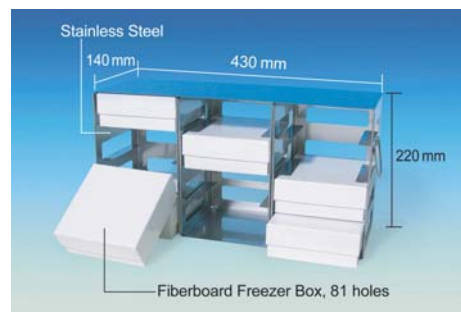

(1) Side Access-type Rack, 20Places "RFU50SA"



(2) Hard Board Box, 81 holes, w/cover



(3) Temp. Chart Recorder, "KT622"



(4) Chart Paper, "C657", 60 charts for 420 days



(5) Replacement Pens 6ea, Red color, "P222", for Cart Recorder

>> **DAIHAN WiRe™ App & Web Service**



7" Full Touch Screen TFT LCD (Smart-Lab™ Controller, Only U80 model)

It is possible to monitor and control the Smart-Lab™ equipment through all smart phone, tablet and web without restraint of time and place.



WiRe™ App & Web Service



(6) LCO2 back up system "LCO2B"



(7) Remote Alarm System "RWAS"

※ See page 6 ~ 11 for the Feature of Smart-Lab™ System